


Computer Controlled **Refrigeration and Air Conditioning Unit**, with SCADA

(two condensers (water and air) and one evaporator (water)



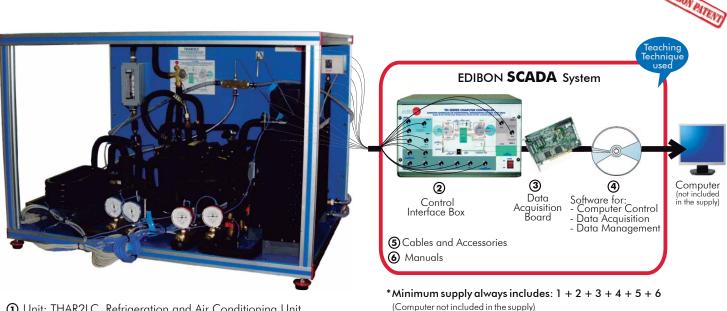

① Unit: THAR2LC. Refrigeration and Air Conditioning Unit (two condensers (water and air) and one evaporator (water))

Key features:

- Advanced Real-Time SCADA.
- Open Control + Multicontrol + Real-Time Control.
- Specialized EDIBON Control Software based on Labview.
- National Instruments Data Acquisition board (250 KS/s , kilo samples per second).
- Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.
- Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.
- Capable of doing applied research, real industrial simulation, training courses, etc.
- Remote operation and control by the user and remote control for EDIBON technical support, are always included.
- Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).
- Designed and manufactured under several quality standards.
- Optional CAL software helps the user perform calculations and comprehend the results.
- This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.

For more information about Key Features, click here:





(total safety



Page 1

You




OPEN CONTROL MULTICONTROL REAL TIME CONTROL





Worlddidac Quality Charter Certificate (Worlddidac Member)

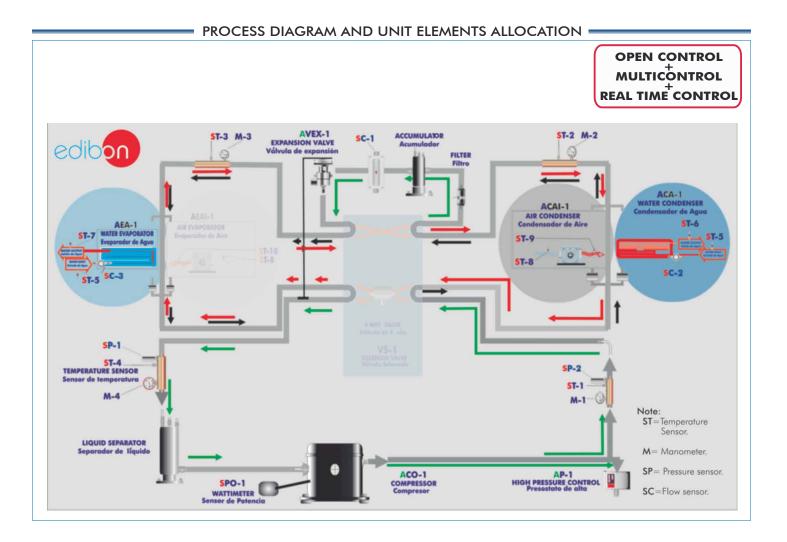

SPECIFICATIONS =

Items supplied as standard

(1) THAR2LC. Unit: Bench-top unit.

Anodized aluminium structure and panels in painted steel.

Main metallic elements in stainless steel.

Diagram in the front panel with similar distribution to the elements in the real unit.

Cooling compressor, computer controlled.

Water condenser. Air condenser, computer controlled.

Water evaporator.

High pressure control.

Coolant accumulation tank.

Cooling filter.

Expansion valve.

Tank of division of the cooling liquid.

4 Manometers.

9 Temperature sensors:

Temperature sensor, J type (compressor outlet).

Temperature sensor, J type (condenser outlet).

Temperature sensor, J type (evaporator inlet).

Temperature sensor, J type (compressor inlet).

Temperature sensor, J type (water inlet).

Temperature sensor, J type (water condenser outlet/water).

Temperature sensor, J type (water evaporator outlet/water).

Temperature sensor, J type (air condenser outlet/air).

Temperature sensor, J type (room air).

3 Flow sensors:

Coolant flow sensor.

Water flow sensor (water condenser). Water flow sensor (water evaporator).

2 Pressure sensors:

Cooling pressure sensor (compressor outlet).

Cooling pressure sensor (compressor inlet).

Wattmeter.

Enthalpy diagram of the refrigerant R134a.

② THAR2LC/CIB. Control Interface Box:

Control interface box with process diagram in the front panel and with the same distribution that the different elements located in the unit, for an easy understanding by the student.

All sensors, with their respective signals, are properly manipulated from -10V. to +10V computer output. Sensors connectors in the interface have different pines numbers (from 2 to 16), to avoid connection errors. Single cable between the control interface box and computer.

The unit control elements are permanently computer controlled, without necessity of changes or connections during the whole process test procedure.

Simultaneous visualization in the computer of all parameters involved in the process. Calibration of all sensors involved in the process.

Real time curves representation about system responses. Storage of all the process data and results in a file. Graphic representation, in real time, of all the process/system responses.

All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process.

All the actuators and sensors values and their responses are displayed on only one screen in the computer.

Shield and filtered signals to avoid external interferences.

Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process. Real time computer control for pumps, compressors, resistances, control valves, etc. Open control allowing modifications, at any moment and in real time, of parameters involved in the process

simultaneously. Three safety levels, one mechanical in the unit, other electronic in the control interface and the third one in the control software.

③ DAB. Data Acquisition Board:

PCI Data acquisition board (National Instruments) to be placed in a computer slot. Bus PCI.

Analog input: Channels=16 single-ended or 8 differential. Resolution=16 bits, 1 in 65536. Sampling rate up to: 250 KS/s (Kilo samples per second). Input range (V)=±10V. Data transfers=DMA, interrupts, programmed I/0. Number of DMA channels=6.

Analog output: Channels=2. Resolution=16 bits, 1 in 65536. Maximum output rate up to: 833 KS/s.

Output range(V)=±10V. Data transfers=DMA, interrupts, programmed I/O. Digital Input/Output: Channels=24 inputs/outputs. D0 or DI Sample Clock frequency: 0 to 1 MHz.

Timing: Counter/timers=2. Resolution: Counter/timers: 32 bits.

④ THAR2LC/CCSOF Computer Control + Data Acquisition + Data Management Software:

Compatible with actual Windows operating systems. Graphic and intuitive simulation of the process in screen. Compatible with the industry standards.

Registration and visualization of all process variables in an automatic and simultaneous way.

Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.

Management, processing, comparison and storage of data.

Sampling velocity up to 250,000 data per second.

Calibration system for the sensors involved in the process.

the minimum supply, enabling a normal operation.

It allows the registration of the alarms state and the graphic representation in real time.

Comparative analysis of the obtained data, after the process and modification of the conditions during the process. Open software, allowing to the teacher to modify texts, instructions. Teacher's and student's passwords to

facilitate the teacher's control on the student, and allowing the access to different work levels.

This unit allows the 30 students of the classroom to visualize simultaneously all results and manipulation of the unit, during the process, by using a projector or an electronic whiteboard.

Cables and Accessories, for normal operation.

Manuals: This unit is **supplied with 8 manuals**: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.



THAR2LC/CIB



DAB



THAR2LC/CCSOF

Continue...

*References 1 to 6: THAR2LC + THAR2LC/CIB + DAB + THAR2LC/CCSOF + Cables and Accessories + Manuals are included in

THAR2LC. Unit

Additional and optional items to the standard supply

PLC. Industrial Control using PLC (7 and 8):

⑦ PLC-PI. PLC Module: Circuit diagram in the front panel. Front panel: Digital inputs(X) and Digital outputs (Y) block: 16 Digital inputs, activated by switches and 16 LEDs for confirmation (red). 14 Digital outputs (through SCSI connector) with 14 LEDs for message (green). Analog inputs block: 16 Analog inputs (-10V. to + 10V.) (through SCSI connector). Analog outputs block: 4 Analog outputs (-10V. to + 10V.) (through SCSI connector). Touch screen: High visibility and multiple functions. Display of a highly visible status. Recipe function. Bar graph function. Flow display function. Alarm list. Multi language function. True type fonts. Back panel: Power supply connector. Fuse 2A. RS-232 connector to PC. USB 2.0 connector to PC. Inside: Power supply outputs: 24 Vdc, 12 Vdc, -12 Vdc, 12 Vdc variable. Panasonic PLC: High-speed scan of 0.32 μ sec. for a basic instruction. Program capacity of 32 Ksteps, with a sufficient comment area. Power supply input (100 to 240 VAC). DC input: 16 (24 V DC). Relay output: 14. High-speed counter. Multi-point PID control. Digital inputs/outputs and analog inputs/outputs Panasonic modules. Communication RS232 wire, to computer (PC). **⑧ THAR2LC/PLC-SOF. PLC Control Software:** For this particular unit, always included with PLC supply.



PLC-PI

Items available on request

THAR2LC/CAL. Computer Aided Learning Software (Results Calculation and Analysis).

⁽ⁱ⁾ THAR2LC/FSS. Faults Simulation System.

Some Practical Possibilities of the Unit:

- 1.-Determination of the inlet power, heat produced and performance coefficient. Water as heat source.
- 2.-Preparation of performance curves of the unit with different inlet and outlet temperatures. Water as heat source.
- 3.-Lay out of the steam compression cycle in a diagram P-H and comparison with the ideal cycle. Water as heat source.
- 4.-Preparation of the performance curves of the unit based on the properties of the refrigerant and at different condensation and evaporation temperatures. Water as heat source.


Other possible practices:

- 5.-Temperature sensors calibration.
- 6.-Flow sensors calibration.
- 7.-Pressure sensors calibration.

Practices to be done by PLC Module (PLC-PI) + PLC Control Software:

- 8.-Control of the THAR2LC unit process through the control interface box without computer.
- 9.-Visualization of all the sensors values used in THAR2LC unit process.
- 10.-Calibration of all sensors included in THAR2LC unit process.
- 11.-Hand on of all the actuators involved in the THAR2LC unit process.
- 12.-Realization of different experiments, in automatic way, without having in front the unit. (This experiment can be decided previously).
- 13.-Simulation of outside actions, in the cases do not exist hardware elements (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).

- 14.-PLC hardware general use and manipulation.
- 15.-PLC process application for the THAR2LC unit.
- 16.-PLC structure.
- 17.-PLC inputs and outputs configuration.
- 18.-PLC configuration possibilities.
- 19.-PLC program languages.
- 20.-PLC different programming standard languages.
- 21.-New configuration and development of new process.
- 22.-Hand on an established process.
- 23.-To visualize and see the results and to make comparisons with the THAR2LC unit process.
- 24.-Possibility of creating new process in relation with the THAR2LC unit.
- 25.-PLC Programming Exercises.
- 26.-Own PLC applications in accordance with teacher and student requirements.

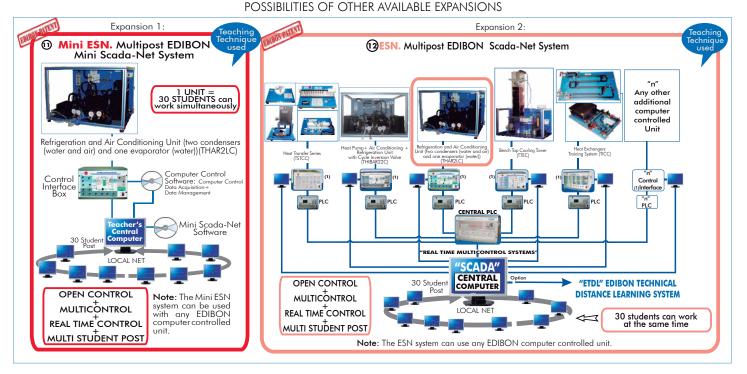

ORDER INFORMATION -

Items supplied as standard:

Minimum configuration for normal operation includes:

- 1 Unit: THAR2LC. Refrigeration and Air Conditioning Unit (two
- condensers (water and air) and one evaporator (water)). ② THAR2LC/CIB. Control Interface Box.
- ③ DAB. Data Acquisition Board.
- THAR2LC/CCSOF. Computer Control + Data Acquisition + Data Management Software.
- **5** Cables and Accessories, for normal operation.
- ⑥ Manuals.

* <u>IMPORTANT:</u> Under <u>THAR2LC</u> we always supply all the elements for immediate running as 1, 2, 3, 4, 5 and 6.

Additional and optional items to the standard supply:

PLC. Industrial Control using PLC (7 and 8):

PCL-PI.PLC Module.

- 8THAR2LC/PLC-SOF. PLC Control Software.
- THAR2LC/CAL. Computer Aided Learning Software (Results Calculation and Analysis). (Available on request).
- OTHAR2LC/FSS. Faults Simulation System. (Available on request).

<u>Expansions</u>

OMini ESN. Multipost EDIBON Mini Scada-Net System.

ESN. Multipost EDIBON Scada-Net System.

REQUIRED SERVICES -

Electrical supply: 220V., 1-phase + neutral + ground, 50 Hz.; or 110V., 1-phase + neutral + ground, 60Hz.; and 1 CV max.

Water supply and drainage.

Computer (PC).

DIMENSIONS & WEIGHTS =

THAR2LC Unit: -Dimensions: 900 x 600 x 500 mm. approx. -Weight: 85 Kg. approx.

Control Interface Box: -Dimensions: 490 x 330 x 310 mm. approx.

-Weight: 10 Kg. approx.

 $\label{eq:PLC-PI} PLC \ Module \ (PLC-PI): \quad \ \ -Dimensions: \ 490 \ x \ 330 \ x \ 310 \ mm. \ approx.$

-Weight: 30 Kg. approx.

RECOMMENDED ACCESSORIES

For refilling R134a refrigerant and maintenance, we recommend:

- T/KIT1. Maintenance Kit, containing: vacuum pump, hoses and manometers.

- T/KIT2. Maintenance Kit, containing: leakage detector.

- R134a refrigerant (to be acquired by the customer locally).

AVAILABLE VERSIONS

Offered in this catalogue:

- THAR2LC. Computer Controlled Refrigeration and Air Conditioning Unit (two condensers (water and air) and one evaporator (water)).

Offered in other catalogue:

- THAR2LB. Refrigeration and Air Conditioning Unit (two condensers (water and air) and one evaporator (water)).

OTHER AVAILABLE REFRIGERATION AND AIR CONDITIONING UNITS -

- THAR22C.	Computer Controlled Refrigeration and Air Conditioning Unit (two condensers (water and air) and two evaporators (water and air)).
- THAR22B.	Refrigeration and Air Conditioning Unit (two condensers (water and air) and two evaporators (water and air)).
- THARL2C.	Computer Controlled Refrigeration and Air Conditioning Unit (one condenser (water) and two evaporators (water and air)).
- THARL2B.	Refrigeration and Air Conditioning Unit (one condenser (water) and two evaporators (water and air)).
- THARALC.	Computer Controlled Refrigeration and Air Conditioning Unit (one condenser (air) and one evaporator (water)).
- THARALB.	Refrigeration and Air Conditioning Unit (one condenser (air) and one evaporator (water)).
- THARA2C.	Computer Controlled Refrigeration and Air Conditioning Unit (one condenser (air) and two evaporators (water and air)).
- THARA2B.	Refrigeration and Air Conditioning Unit (one condenser (air) and two evaporators (water and air)).
- THARLLC.	Computer Controlled Refrigeration and Air Conditioning Unit (one condenser (water) and one evaporator (water)).
- THARLLB.	Refrigeration and Air Conditioning Unit (one condenser (water) and one evaporator (water)).

*Specifications subject to change without previous notice, due to the convenience of improvements of the product.



C/ Del Agua, 14. Polígono Industrial San José de Valderas. 28918 LEGANÉS. (Madrid). SPAIN. Phone: 34-91-6199363 FAX: 34-91-6198647 E-mail: edibon@edibon.com WEB site: **www.edibon.com**

Issue: ED01/13 Date: May/2013 REPRESENTATIVE: