

* Minimum supply always includes: 1 + 2 + 3 + 4 + 5 + 6
(Computer not included in the supply)

① Unit: TCEC. Separating & Throttling Calorimeter

Key features:

- ▶ **Advanced Real-Time SCADA.**
- ▶ **Open Control + Multicontrol + Real-Time Control.**
- ▶ **Specialized EDIBON Control Software based on Labview.**
- ▶ **National Instruments Data Acquisition board (250 KS/s , kilo samples per second).**
- ▶ **Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.**
- ▶ **Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.**
- ▶ **Capable of doing applied research, real industrial simulation, training courses, etc.**
- ▶ **Remote operation and control by the user and remote control for EDIBON technical support, are always included.**
- ▶ **Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).**
- ▶ **Designed and manufactured under several quality standards.**
- ▶ **Optional CAL software helps the user perform calculations and comprehend the results.**
- ▶ **This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.**

**OPEN CONTROL
+
MULTICONTROL
+
REAL TIME CONTROL**

For more information about Key Features, click here:



ISO 9000: Quality Management
(for Design, Manufacturing,
Commercialization and After-sales service)



European Union Certificate
(total safety)



Certificates ISO 14000 and
ECO-Management and Audit Scheme
(environmental management)



Worlddidac Quality Charter
Certificate
(Worlddidac Member)

www.edibon.com

- ↳ Products
- ↳ Products range
- ↳ Units
- ↳ 9 - Thermodynamics & Thermotechnics

GENERAL DESCRIPTION

Unit, computer controlled, to determine low and high water contents in two-phase liquid-water mixture and the dryness fraction of steam by means a separating and throttling calorimeters.


TCESC unit makes it possible to determine the content of water in the wet steam, both in big and small quantities. Such steam is generated in a steam generator or in a boiler.

The separating calorimeter is used for a direct determination of great quantities of water (up to 0.5). In case of small contents of humidity (the quality of the steam ranges from 0.95 to 1.0) can be determined indirectly with the throttling calorimeter, by means of the adiabatic expansion and the conversion into superheated steam, with the aid of the h-s diagram.

At the same time, the TCESC unit includes two heat exchangers for each condensate obtained in both calorimeters. The condenser belonging to the throttling calorimeter makes possible the study of the heat transfer between the condensate of such calorimeter (hot water) which flows through a coil and the cold water that flows in the exchanger. Such condenser makes possible the measurement of the temperature at the inlet and outlet of the exchanger, both in the cold and hot water.

This Computer Controlled Unit is supplied with the EDIBON Computer Control System (SCADA), and includes: The unit itself + a Control Interface Box + a Data Acquisition Board + Computer Control and Data Acquisition Software Packages, for controlling the process and all parameters involved in the process.

PROCESS DIAGRAM AND UNIT ELEMENTS ALLOCATION

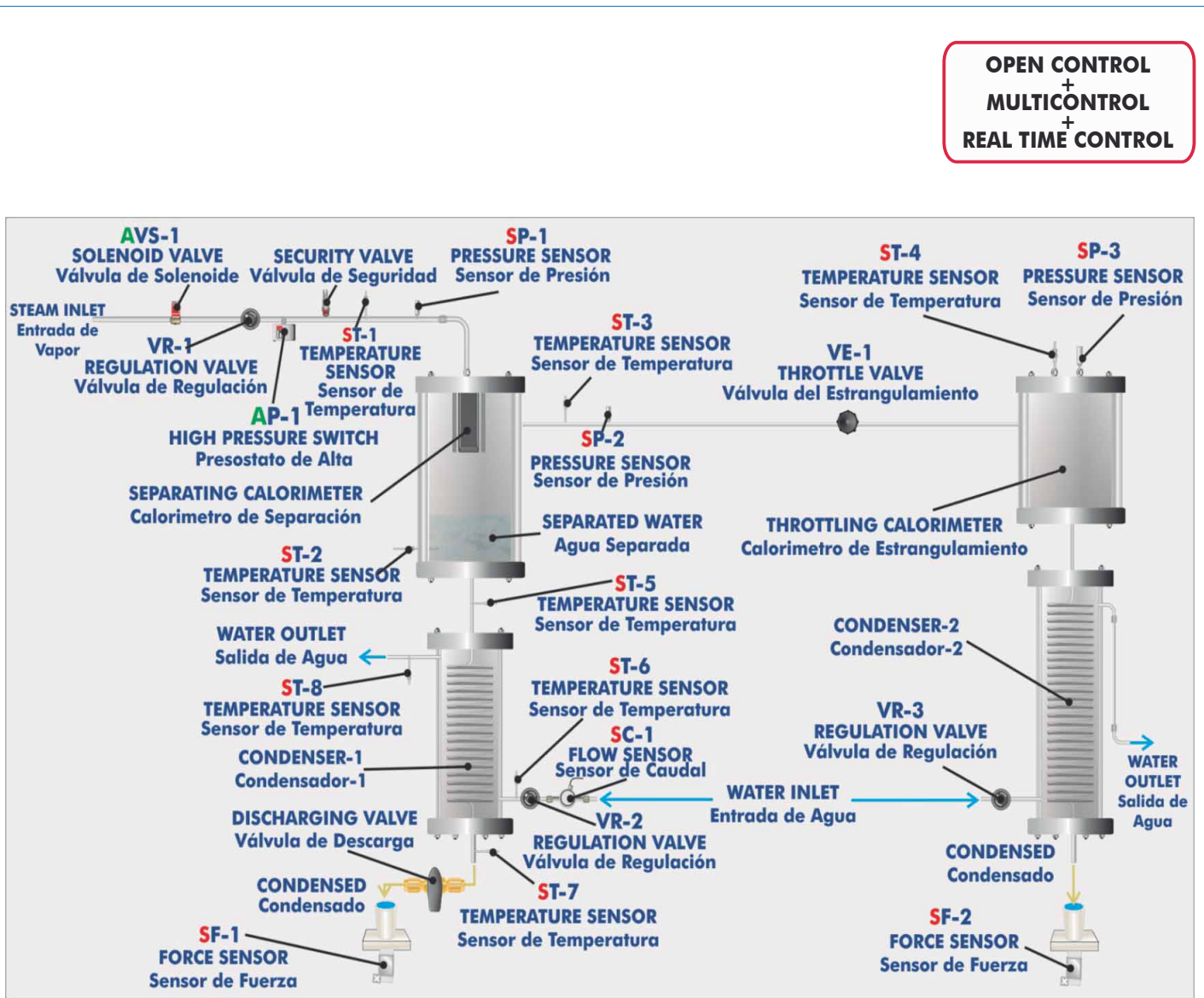

With this unit there are several options and possibilities:

- Main items: 1, 2, 3, 4, 5 and 6.
- Optional items: 7, 8, 9, 10, 11 and 12.

Let us describe first the main items (1 to 6):

① TCESC. Unit:

Anodized aluminium structure and panels in painted steel.

Main metallic elements in stainless steel.

Diagram in the front panel with similar distribution to the elements in the real unit.

Separating calorimeter, with a water level sight glass, for a direct determination of big contents of water in two-phase liquid-water mixture.

Throttling calorimeter for the determination of small humidity contents.

Water condenser connected to the separating calorimeter. It consists of a stainless steel vessel and a copper coil:

Vessel inner diameter: 70 mm.

Vessel length: 160 mm.

Coil inner diameter: 4.35 mm.

Coil external diameter: 6.35 mm.

A water condenser connected to the throttling calorimeter. It consists of a stainless steel vessel and a copper coil:

Vessel inner diameter: 80 mm.

Vessel length: 360 mm.

Coil inner diameter: 4.35 mm.

Coil external diameter: 6.35 mm.

8 "J type" temperature sensors:

4 Temperature sensors to determine the steam quality.

2 Temperature sensors to measure the temperature of the hot water (coming from the throttling calorimeter).

2 Temperature sensors to measure the temperature of the cold water.

3 Pressure sensors, range: 0 -10 bar.

Water flow sensor, range: 0.25 - 6.5 l./min.

Two force sensors to measure the weight of condensed water in both calorimeters. Range of the sensors: 0 - 0.6 Kg. and 0 - 2 Kg.

Two graduated glass vessels, capacity: 250 ml and 2 liters.

High pressure switch.

Safety valve.

2 Solenoid valves.

The complete unit includes as well:

Advanced Real-Time SCADA.

Open Control + Multicontrol + Real-Time Control.

Specialized EDIBON Control Software based on Labview.

National Instruments Data Acquisition board (250 KS/s , kilo samples per second).

Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.

Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.

Capable of doing applied research, real industrial simulation, training courses, etc.

Remote operation and control by the user and remote control for EDIBON technical support, are always included.

Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).

Designed and manufactured under several quality standards.

Optional CAL software helps the user perform calculations and comprehend the results.

This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.



TCESC. Unit

② **TCESC/CIB. Control Interface Box:**

The Control Interface Box is part of the SCADA system.

Control interface box with process diagram in the front panel and with the same distribution that the different elements located in the unit, for an easy understanding by the student.

All sensors, with their respective signals, are properly manipulated from -10V. to +10V. computer output. Sensors connectors in the interface have different pines numbers (from 2 to 16), to avoid connection errors. Single cable between the control interface box and computer.

The unit control elements are permanently computer controlled, without necessity of changes or connections during the whole process test procedure.

Simultaneous visualization in the computer of all parameters involved in the process.

Calibration of all sensors involved in the process.

Real time curves representation about system responses.

Storage of all the process data and results in a file.

Graphic representation, in real time, of all the process/system responses.

All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process. All the actuators and sensors values and their responses are displayed on only one screen in the computer.

Shield and filtered signals to avoid external interferences.

Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process. Real time computer control for pumps, compressors, resistances, control valves, etc. Real time computer control for parameters involved in the process simultaneously.

Open control allowing modifications, at any moment and in real time, of parameters involved in the process simultaneously.

Three safety levels, one mechanical in the unit, another electronic in the control interface and the third one in the control software.



TCESC/CIB

③ **DAB. Data Acquisition Board:**

The Data Acquisition board is part of the SCADA system.

PCI Data acquisition board (National Instruments) to be placed in a computer slot. Bus PCI.

Analog input: Number of channels= 16 single-ended or 8 differential. Resolution= 16 bits, 1 in 65536.

Sampling rate up to: 250 KS/s (kilo samples per second). Input range (V) = ±10 V.

Data transfers=DMA, interrupts, programmed I/O. DMA channels=6.

Analog output: Number of channels=2. Resolution= 16 bits, 1 in 65536.

Maximum output rate up to: 833 KS/s.

Output range(V) = ±10 V. Data transfers=DMA, interrupts, programmed I/O.

Digital Input/Output: Number of channels=24 inputs/outputs.

D0 or DI Sample Clock frequency: 0 to 1 MHz.

Timing: Number of Counter/timers=2. Resolution: Counter/timers: 32 bits.



DAB

④ **TCESC/CCSOF. Computer Control + Data Acquisition + Data Management Software:**

The three softwares are part of the SCADA system.

Compatible with actual Windows operating systems. Graphic and intuitive simulation of the process in screen. **Compatible with the industry standards.**

Registration and visualization of all process variables in an automatic and simultaneous way.

Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.

Management, processing, comparison and storage of data.

Sampling velocity up to 250 KS/s (kilo samples per second).


Calibration system for the sensors involved in the process.

It allows the registration of the alarms state and the graphic representation in real time.

Comparative analysis of the obtained data, after the process and modification of the conditions during the process.

Open software, allowing the teacher to modify texts, instructions. Teacher's and student's passwords to facilitate the teacher's control on the student, and allowing the access to different work levels.

This unit allows the 30 students of the classroom to visualize simultaneously all the results and the manipulation of the unit, during the process, by using a projector or an electronic whiteboard.

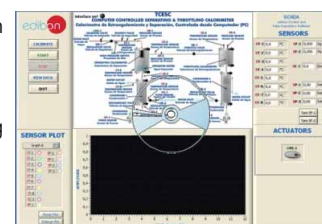

TCESC/CCSOF

⑤ **Cables and Accessories**, for normal operation.

⑥ **Manuals:** This unit is supplied with 8 manuals: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.

* References 1 to 6 are the main items: TCESC + TCESC/CIB + DAB + TCESC/CCSOF + Cables and Accessories + Manuals are included in the minimum supply for enabling normal and full operation.

EXERCISES AND PRACTICAL POSSIBILITIES TO BE DONE WITH MAIN ITEMS

- 1.- To use a separating calorimeter for high water contents.
- 2.- To use a throttling calorimeter for high vapour contents.
- 3.- Determination of the dryness fraction of steam.
- 4.- Global energy balance in the exchanger and study of losses.
- 5.- Determination of the exchanger efficiency. NTU method.

Additional practical possibilities:

- 6.- Sensors calibration.

Other possibilities to be done with this Unit:

- 7.- Many students view results simultaneously.

To view all results in real time in the classroom by means of a projector or an electronic whiteboard.

- 8.- Open Control, Multicontrol and Real Time Control.

This unit allows intrinsically and/or extrinsically to change the span, gains; proportional, integral, derivate parameters; etc, in real time.

- 9.- The Computer Control System with SCADA allows a real industrial simulation.

- 10.- This unit is totally safe as uses mechanical, electrical and electronic, and software safety devices.

- 11.- This unit can be used for doing applied research.

- 12.- This unit can be used for giving training courses to Industries even to other Technical Education Institutions.

- 13.- Control of the TCESC unit process through the control interface box without the computer.

- 14.- Visualization of all the sensors values used in the TCESC unit process.

- By using PLC-PI additional 19 more exercises can be done.

- Several other exercises can be done and designed by the user.

REQUIRED SERVICES

- Electrical supply: single-phase, 220 V./50 Hz. or 110 V./60 Hz.
- Computer (PC).
- Water supply for the service lines to the condensers.
- Water drain for the service lines to the condensers.
- Steam supply, with a maximum working pressure of 10 bar and a temperature of 240°C. EDIBON can offer to work with this unit: TVG. Steam Generator (3 kW) or TGV-6kW. Steam Generator (6kW).
- Chronometer.

DIMENSIONS & WEIGHTS

TCESC:

Unit: -Dimensions: 1000 x 870 x 1590 mm. approx.
(39.37 x 34.25 x 62.59 inches approx.)

-Weight: 70 Kg. approx.
(154.3 pounds approx.)

Control Interface Box:-Dimensions: 490 x 330 x 310 mm. approx.
(19.29 x 12.99 x 12.2 inches approx.)

-Weight: 10 Kg. approx.
(22 pounds approx.)

SCADA
Main screen

The SCADA main screen is divided into several functional areas:

- Top Left:** Edibon logo and a vertical menu with buttons: CALIBRATE, START, STOP, VIEW DATA, QUIT.
- Top Center:** A detailed process diagram of a 'COMPUTER CONTROLLED SEPARATING & THROTTLING CALORIMETER'. It includes components like Steam Inlet, Solenoid Valve (AVS-1), Regulation Valve (VR-1), High Pressure Switch, Separating Calorimeter, Throttling Calorimeter, Condenser-2, Water Inlet, and Water Outlet. Various sensors (ST-1 to ST-8, SF-1, SF-2, SC-1, SP-1, SP-2, SP-3) and valves (VE-1, VR-2, VR-3) are indicated.
- Top Right:** 'SENSORS' display showing real-time values for 11 different sensors (ST-1 to ST-8, SF-1 to SF-2, SC-1, SP-1 to SP-3) with their respective units (°C, kg, l/min, bar).
- Bottom Left:** 'SENSOR PLOT' section with a 'Graph A' dropdown and a list of sensors to be plotted (ST-1 to ST-8, SC-1, SF-1).
- Bottom Center:** A real-time graph showing 'AMPLITUDE' vs 'TIME (seconds)'. The plot area is currently empty.
- Bottom Right:** 'ACTUATORS' section showing a control for the AVS-1 solenoid valve.

- Ⓘ Main software operation possibilities.
- Ⓜ Sensors displays, real time values, and extra output parameters. Sensors: ST=Temperature sensor. SF=Force sensor. SC=Flow sensor. SP=Pressure sensor.
- Ⓝ Actuators controls. Actuator: AVS=Solenoid valve.
- Ⓥ Channel selection and other plot parameters.
- Ⓟ Real time graphics displays.

Software for Sensors Calibration

The calibration software consists of two main windows:

- CALIBRATION:** A window for setting up a single sensor. It shows 'Analog Input Channel' (ST-6), 'Sensor Name' (ST-6), 'Calibration units' (°C), 'Full Scale' (700), 'Gain' (99,1955), and 'Offset' (0,29761). It includes a 'Least Squares Fit' button and 'ENTER', 'EXIT', and 'EXIT & SAVE' buttons.
- MULTICALIBRATE:** A window for calibrating multiple sensors. It features a 'Reference Value' (23,2113), 'AT Full Scale' (100), and 'Tolerance (%)' (1). A table lists sensors and their calibration data:

Reference Select	Sensors	Volts	Calibrated	Err (%)
<input checked="" type="checkbox"/>	ST-1	0,2046	22,3821	0,82
<input checked="" type="checkbox"/>	ST-2	0,2292	23,483	0,28
<input checked="" type="checkbox"/>	ST-3	0,2353	23,1522	0,05
<input checked="" type="checkbox"/>	ST-4	0,2301	23,2113	0,01
<input type="checkbox"/>		0,1527	13,1629	10,04
<input type="checkbox"/>	SEC-3	-5,2792	172,5164	149,31
<input type="checkbox"/>		-0,2362	-22,6609	45,87
<input type="checkbox"/>	SC-1	-0,1774	0,0319629	23,17
<input type="checkbox"/>		-0,2681	-40,4623	83,67
<input type="checkbox"/>		-0,2251	0,4208	22,79
<input type="checkbox"/>		-0,2529	-0,2529	23,46
<input type="checkbox"/>		-0,2063	-0,1178	23,32
<input type="checkbox"/>		-0,2581	-226,9384	250,14
<input type="checkbox"/>		-0,3634	-0,3634	23,57
<input type="checkbox"/>		-0,275	-0,275	23,48
<input type="checkbox"/>		-0,2005	-0,2005	23,41


 The window also includes 'GAIN', 'OFFSET', and 'p' columns for each sensor and 'ENTER', 'DONE' buttons.

By using a free of charge code, the teacher and the students can calibrate the unit.

The teacher can recover his/her own calibration by using the EDIBON code that we give free of charge.

SOME TYPICAL RESULTS

We see the sensors values and graphic representation in real time.



START


STOP

SAVING DATA

Periodic (seconds)
1,9
6,2 sec

Take Data

TCESC
COMPUTER CONTROLLED SEPARATING & THROTTLING CALORIMETER
Calorímetro de Estrangulamiento y Separación, Controlada desde Computador (PC)

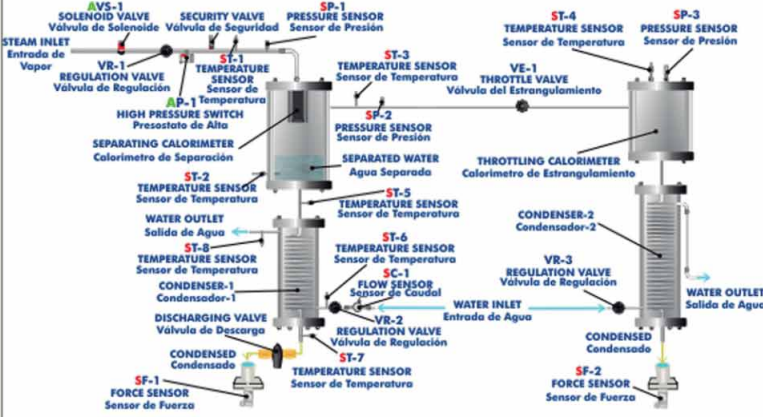

SCADA
edibon Control and
Data Acquisition Software

SENSORS

ST-1	24,4 °C	SF-1	0,000 kg
ST-2	26,2 °C	SF-2	-0,705 kg
ST-3	26,1 °C	SC-1	0,0 l/min
ST-4	26,9 °C	SP-1	0,00 bar
ST-5	24,7 °C	SP-2	0,00 bar
ST-6	23,7 °C	SP-3	0,00 bar
ST-7	0,0 °C		
ST-8	0,3 °C		

Tare SF-1
Tare SF-2

ACTUATORS

AVS-1


SENSOR PLOT

Graph A

- ST-1 SF-2
- ST-2 SP-1
- ST-3 SP-2
- ST-4 SP-3
- ST-5
- ST-6
- ST-7
- ST-8
- SC-1
- SF-1

Reset Plot
Enlarge Plot


Graph A

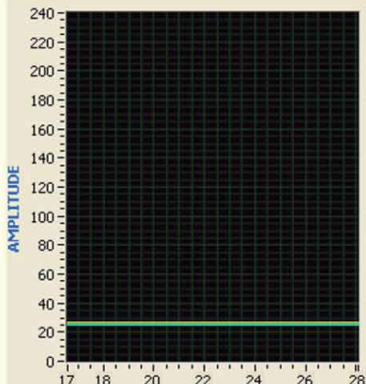

AMPLITUDE

TIME (seconds)

Graph B

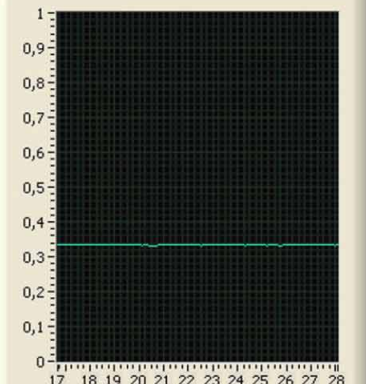

AMPLITUDE

TIME (seconds)


VIEWDATA=Graphic visualization of the saved data in a file.

VIEWDATA

View Data Select

<input checked="" type="checkbox"/>	ST-1	<input type="checkbox"/>	1	▽	ST-1
<input checked="" type="checkbox"/>	ST-2	<input type="checkbox"/>	1	▽	ST-2
<input checked="" type="checkbox"/>	ST-3	<input type="checkbox"/>	1	▽	ST-3
<input checked="" type="checkbox"/>	ST-4	<input type="checkbox"/>	1	▽	ST-4
<input checked="" type="checkbox"/>	ST-5	<input type="checkbox"/>	1	▽	ST-5
<input checked="" type="checkbox"/>	ST-6	<input type="checkbox"/>	1	▽	ST-6
<input checked="" type="checkbox"/>	ST-7	<input type="checkbox"/>	1	▽	ST-7
<input checked="" type="checkbox"/>	ST-8	<input type="checkbox"/>	1	▽	ST-8
<input type="checkbox"/>	SC-1	<input type="checkbox"/>	1	▽	SC-1
<input type="checkbox"/>	SF-1	<input type="checkbox"/>	1	▽	SF-1
<input type="checkbox"/>	SF-2	<input type="checkbox"/>	1	▽	SF-2
<input checked="" type="checkbox"/>	SP-1	<input type="checkbox"/>	1	▽	SP-1
<input type="checkbox"/>	SP-2	<input type="checkbox"/>	1	▽	SP-2
<input type="checkbox"/>	SP-3	<input type="checkbox"/>	1	▽	SP-3
<input type="checkbox"/>	AVS-1	<input type="checkbox"/>	1	▽	AVS-1

Data Graph

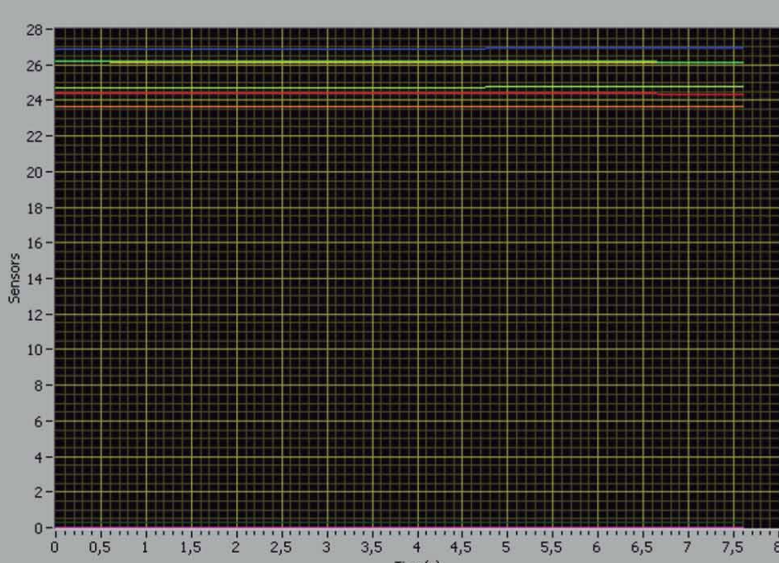

Sensors

Time(s)

AutoScale

CLOSE

COMPLETE TECHNICAL SPECIFICATIONS (for optional items)

Additionally to the main items (1 to 6) described, we can offer, as optional, other items from 7 to 12.

All these items try to give more possibilities for:

- a) Industrial configuration. (PLC)
- b) Technical and Vocational Education configuration. (CAL and FSS)
- c) Higher Education and/or Technical and Vocational Education configuration. (CAL)
- d) Multipost Expansions options. (Mini ESN and ESN)

a) Industrial configuration

⑦ **PLC. Industrial Control using PLC** (it includes PLC-PI Module plus PLC-SOF Control Software):

-PLC-PI. PLC Module:

Metallic box.

Circuit diagram in the module front panel.

Front panel:

Digital inputs(X) and Digital outputs (Y) block:

16 Digital inputs, activated by switches and 16 LEDs for confirmation (red).

14 Digital outputs (through SCSI connector) with 14 LEDs for message (green).

Analog inputs block:

16 Analog inputs (- 10 V. to + 10 V.) (through SCSI connector).

Analog outputs block:

4 Analog outputs (- 10 V. to + 10 V.) (through SCSI connector).

Touch screen:

High visibility and multiple functions. Display of a highly visible status. Recipe function. Bar graph function. Flow display function. Alarm list.

Multi language function. True type fonts.

Back panel:

Power supply connector. Fuse 2A. RS-232 connector to PC. USB 2.0 connector to PC.

Inside:

Power supply outputs: 24 Vdc, 12 Vdc, -12 Vdc, 12 Vdc variable.

Panasonic PLC:

High-speed scan of 0.32 μsec. for a basic instruction.

Program capacity of 32 Ksteps, with a sufficient comment area.

Power supply input (100 to 240 V AC).

DC input: 16 (24 V DC).

Relay output: 14.

High-speed counter.

Multi-point PID control.

Digital inputs/outputs and analog inputs/outputs Panasonic modules.


Communication RS232 wire to computer (PC).

Dimensions: 490 x 330 x 310 mm. approx. (19.29 x 12.99 x 12.2 inches approx.). Weight: 30 Kg. approx. (66 pounds approx.).

-TCESC/PLC-SOF. PLC Control Software:

For this particular unit, always included with PLC supply.

The software has been designed using Labview and it follows the unit operation procedure and linked with the Control Interface Box used in the Computer Controlled Separating & Throttling Calorimeter (TCESC).

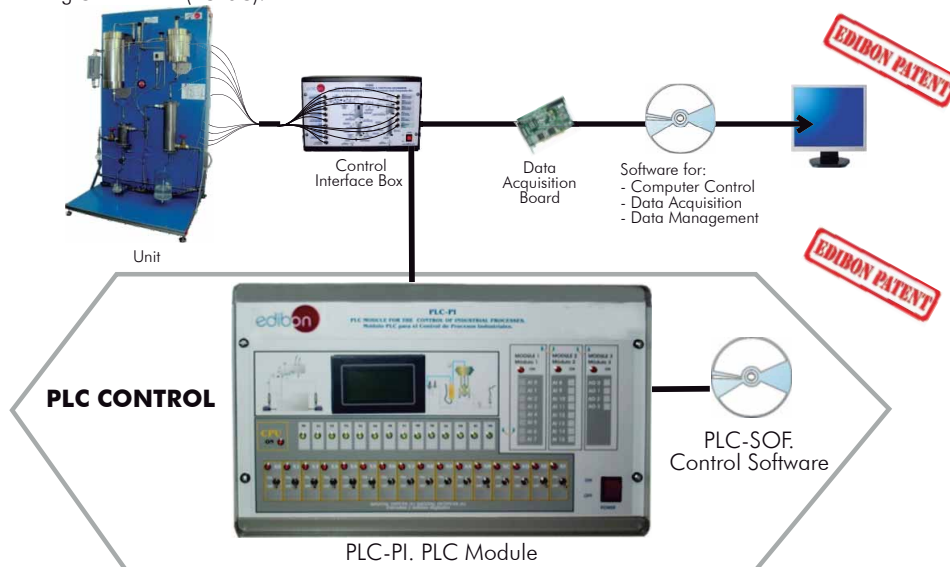

Practices to be done with PLC-PI:

- 1.- Control of the TCESC unit process through the control interface box without the computer.
- 2.- Visualization of all the sensors values used in the TCESC unit process.
- 3.- Calibration of all sensors included in the TCESC unit process.
- 4.- Hand on of all the actuators involved in the TCESC unit process.
- 5.- Realization of different experiments, in automatic way, without having in front the unit. (This experiment can be decided previously).
- 6.- Simulation of outside actions, in the cases hardware elements do not exist. (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).
- 7.- PLC hardware general use and manipulation.
- 8.- PLC process application for TCESC unit.
- 9.- PLC structure.
- 10.- PLC inputs and outputs configuration.
- 11.- PLC configuration possibilities.
- 12.- PLC programming languages.
- 13.- PLC different programming standard languages.
- 14.- New configuration and development of new process.
- 15.- Hand on an established process.
- 16.- To visualize and see the results and to make comparisons with the TCESC unit process.
- 17.- Possibility of creating new process in relation with the TCESC unit.
- 18.- PLC Programming exercises.
- 19.- Own PLC applications in accordance with teacher and student requirements.

b) Technical and Vocational Education configuration

⑧ **TCESC/CAI. Computer Aided Instruction Software System.**

This complete package includes two Softwares: the INS/SOF. Classroom Management Software (Instructor Software) and the TCESC/SOF. Computer Aided Instruction Software (Student Software).

This software is optional and can be used additionally to items (1 to 6).

This complete package consists on an Instructor Software (INS/SOF) totally integrated with the Student Software (TCESC/SOF). Both are interconnected so that the teacher knows at any moment what is the theoretical and practical knowledge of the students. These, on the other hand, get a virtual instructor who helps them to deal with all the information on the subject of study.

- INS/SOF. Classroom Management Software (Instructor Software):

The Instructor can:

- Organize Students by Classes and Groups.
- Create easily new entries or delete them.
- Create data bases with student information.
- Analyze results and make statistical comparisons.
- Print reports.
- Develop own examinations.
- Detect student's progress and difficulties.
- ...and many other facilities.

Instructor Software



- TCESC/SOF. Computer Aided Instruction Software (Student Software):

It explains how to use the unit, run the experiments and what to do at any moment.

This Software contains:

- Theory.
- Exercises.
- Guided Practices.
- Exams.

Student Software



For more information see CAI catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/CAI.pdf

⑨ **TCESC/FSS. Faults Simulation System.**

Faults Simulation System (FSS) is a Software package that simulates several faults in any EDIBON Computer Controlled Unit. It is useful for Technical and Vocational level.

The "FAULTS" mode consists on causing several faults in the unit normal operation. The student must find them and solve them.

There are several kinds of faults that can be grouped in the following sections:

Faults affecting the sensors measurement:

- An incorrect calibration is applied to them.
- Non-linearity.

Faults affecting the actuators:

- Actuators channels interchange at any time during the program execution.
- Response reduction of an actuator.


Faults in the controls execution:

- Inversion of the performance in ON/OFF controls.
- Reduction or increase of the calculated total response.
- The action of some controls is annulled.

On/off faults:

- Several on/off faults can be included.

Example of some screens

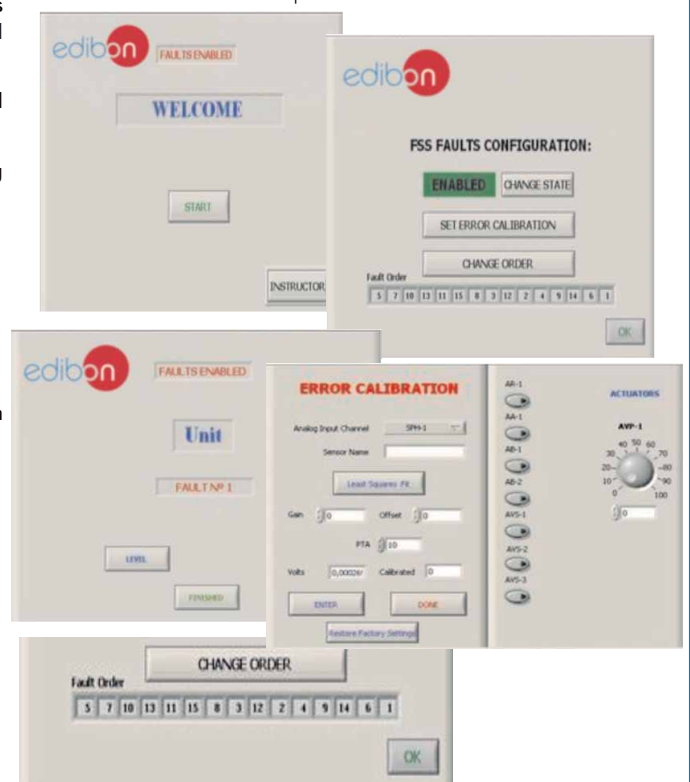

For more information see FSS catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/FSS.pdf

c) Higher Education and/or Technical and Vocational Education configuration

⑩ **TCESC/CAL. Computer Aided Learning Software (Results Calculation and Analysis).**

This Computer Aided Learning Software (CAL) is a Windows based software, simple and very easy to use, specifically developed by EDIBON. It is very useful for Higher Education level.

CAL is a class assistant that helps in making the necessary calculations to extract the right conclusions from data obtained during the experimental practices.

CAL will perform the calculations.

CAL computes the value of all the variables involved.


It allows to plot and print the results. Between the plotting options, any variable can be represented against any other.

Different plotting displays.

It has a wide range of information, such as constant values, unit conversion factors and integral and derivative tables.

For more information see CAL catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/CAL.pdf

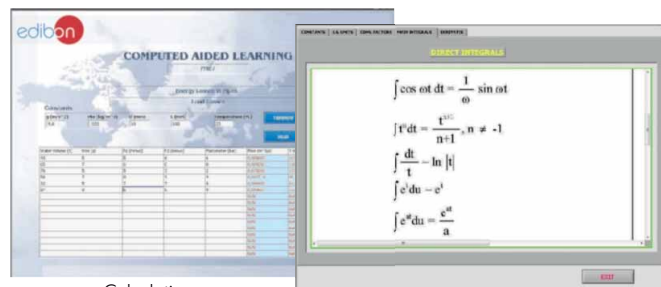

Calculations

Information of constant values, unit conversion factors and integral and derivative tables



Plotting options

d) Multipost Expansions options

⑪ **Mini ESN. EDIBON Mini Scada-Net System.**

Mini ESN. EDIBON Mini Scada-Net System allows up to 30 students to work with a Teaching Unit in any laboratory, simultaneously. It is useful for both, Higher Education and/or Technical and Vocational Education.

The Mini ESN system consists on the adaptation of any EDIBON Computer Controlled Unit with SCADA integrated in a local network.

This system allows to view/control the unit remotely, from any computer integrated in the local net (in the classroom), through the main computer connected to the unit. Then, the number of possible users who can work with the same unit is higher than in an usual way of working (usually only one).

Main characteristics:


- It allows up to 30 students to work simultaneously with the EDIBON Computer Controlled Unit with SCADA, connected in a local net.
- Open Control + Multicontrol + Real Time Control + Multi Student Post.
- Instructor controls and explains to all students at the same time.
- Any user/student can work doing "real time" control/multicontrol and visualisation.
- Instructor can see in the computer what any user/student is doing in the unit.
- Continuous communication between the instructor and all the users/students connected.

Main advantages:

- It allows an easier and quicker understanding.
- This system allows you can save time and cost.
- Future expansions with more EDIBON Units.

For more information see Mini ESN catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/Mini-ESN.pdf

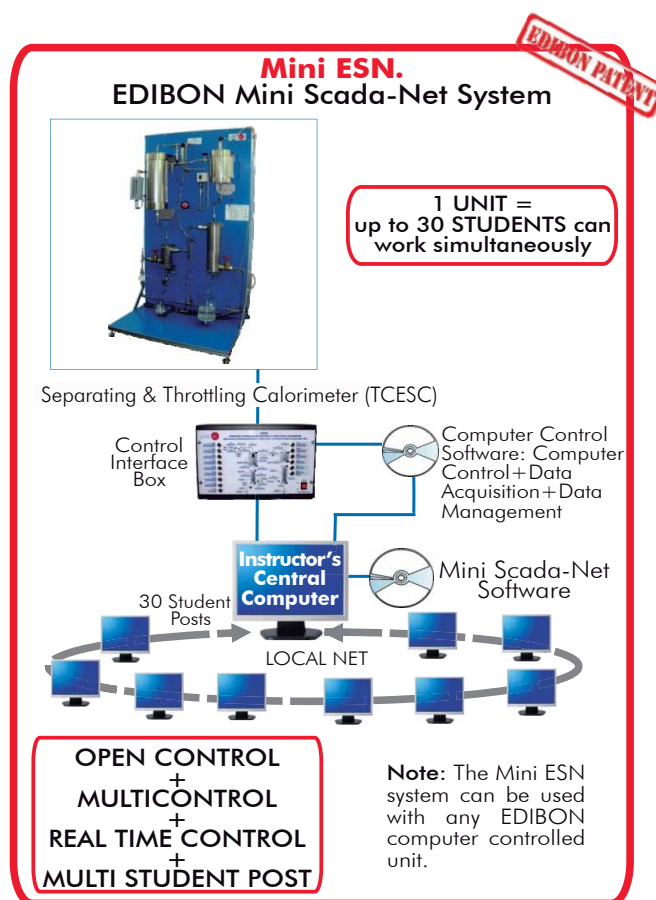

⑫ **ESN. EDIBON Scada-Net System.**

This unit can be integrated, in future, in a Complete Laboratory with many Units and many Students.

For more information see ESN catalogue. Click on the following link:

www.edibon.com/products/catalogues/en/units/thermodynamicsthermotechnics/esn-thermodynamics/ESN-THERMODYNAMICS.pdf

ORDER INFORMATION

Main items (always included in the supply)

Minimum supply always includes:

- ① **Unit: TCESC. Separating & Throttling Calorimeter.**
- ② **TCESC/CIB. Control Interface Box.**
- ③ **DAB. Data Acquisition Board.**
- ④ **TCESC/CCSOF. Computer Control + Data Acquisition + Data Management Software.**
- ⑤ **Cables and Accessories**, for normal operation.
- ⑥ **Manuals.**

* **IMPORTANT:** Under TCESC we always supply all the elements for immediate running as 1, 2, 3, 4, 5 and 6.

Optional items (supplied under specific order)

a) Industrial configuration

- ⑦ PLC. Industrial Control using PLC (it includes PLC-PI Module plus PLC-SOF Control Software):
 - PCL-PI. PLC Module.
 - TCESC/PLC-SOF. PLC Control Software.

b) Technical and Vocational configuration

- ⑧ TCESC/CAI. Computer Aided Instruction Software System.
- ⑨ TCESC/FSS. Faults Simulation System.

c) Higher Education and/or Technical and Vocational Education configuration

- ⑩ TCESC/CAL. Computer Aided Learning Software (Results Calculation and Analysis).

d) Multipost Expansions options

- ⑪ Mini ESN. EDIBON Mini Scada-Net System.
- ⑫ ESN. EDIBON Scada-Net System.

① TCESC. Unit:

Anodized aluminium structure and panels in painted steel.
 Main metallic elements in stainless steel.
 Diagram in the front panel with similar distribution to the elements in the real unit.
 Separating calorimeter, with a water level sight glass, for a direct determination of big contents of water in two-phase liquid-water mixture.
 Throttling calorimeter for the determination of small humidity contents.
 Water condenser connected to the separating calorimeter. It consists of a stainless steel vessel and a copper coil:
 Vessel inner diameter: 70 mm.
 Vessel length: 160 mm.
 Coil inner diameter: 4.35 mm.
 Coil external diameter: 6.35 mm.

A water condenser connected to the throttling calorimeter. It consists of a stainless steel vessel and a copper coil:
 Vessel inner diameter: 80 mm.
 Vessel length: 360 mm.
 Coil inner diameter: 4.35 mm.
 Coil external diameter: 6.35 mm.

8 " J type" temperature sensors:
 4 Temperature sensors to determine the steam quality.
 2 Temperature sensors to measure the temperature of the hot water (coming from the throttling calorimeter).
 2 Temperature sensors to measure the temperature of the cold water.

3 Pressure sensors, range: 0 - 10 bar.
 Water flow sensor, range: 0.25 - 6.5 l./min.
 Two force sensors to measure the weight of condensed water in both calorimeters. Range of the sensors: 0 - 0.6 Kg. and 0 - 2 Kg.
 Two graduated glass vessels, capacity: 250 ml and 2 liters.
 High pressure switch.
 Safety valve.
 2 Solenoid valves.

The complete unit includes as well:
 Advanced Real-Time SCADA.
 Open Control + Multicontrol + Real-Time Control.
 Specialized EDIBON Control Software based on Labview.
 National Instruments Data Acquisition board (250 KS/s , kilo samples per second).
 Calibration exercises, which are included, teach the user how to calibrate a sensor and the importance of checking the accuracy of the sensors before taking measurements.
 Projector and/or electronic whiteboard compatibility allows the unit to be explained and demonstrated to an entire class at one time.
 Capable of doing applied research, real industrial simulation, training courses, etc.
 Remote operation and control by the user and remote control for EDIBON technical support, are always included.
 Totally safe, utilizing 4 safety systems (Mechanical, Electrical, Electronic & Software).
 Designed and manufactured under several quality standards.
 Optional CAL software helps the user perform calculations and comprehend the results.
 This unit has been designed for future expansion and integration. A common expansion is the EDIBON Scada-Net (ESN) System which enables multiple students to simultaneously operate many units in a network.

② TCESC/CIB. Control Interface Box:

The Control Interface Box is part of the SCADA system. Control interface box with process diagram in the front panel.
 The unit control elements are permanently computer controlled.
 Simultaneous visualization in the computer of all parameters involved in the process.
 Calibration of all sensors involved in the process.
 Real time curves representation about system responses.
 All the actuators' values can be changed at any time from the keyboard allowing the analysis about curves and responses of the whole process.
 Shield and filtered signals to avoid external interferences.
 Real time computer control with flexibility of modifications from the computer keyboard of the parameters, at any moment during the process.
 Real time computer control for parameters involved in the process simultaneously.
 Open control allowing modifications, at any moment and in real time, of parameters involved in the process simultaneously.
 Three safety levels, one mechanical in the unit, another electronic in the control interface and the third one in the control software.

③ DAB. Data Acquisition Board:

The Data Acquisition board is part of the SCADA system.
 PCI Data acquisition board (National Instruments) to be placed in a computer slot.
 Analog input: Channels= 16 single-ended or 8 differential. Resolution= 16 bits, 1 in 65536. Sampling rate up to: 250 KS/s (kilo samples per second).
 Analog output: Channels=2. Resolution= 16 bits, 1 in 65536.
 Digital Input/Output: Channels=24 inputs/outputs.

④ TCESC/CCSOF. Computer Control + Data Acquisition + Data Management Software:

The three softwares are part of the SCADA system.
 Compatible with the industry standards.
 Flexible, open and multicontrol software, developed with actual windows graphic systems, acting simultaneously on all process parameters.
 Management, processing, comparison and storage of data.
 Sampling velocity up to 250 KS/s (kilo samples per second).
 Calibration system for the sensors involved in the process.
 It allows the registration of the alarms state and the graphic representation in real time.
 Open software, allowing the teacher to modify texts, instructions. Teacher's and student's passwords to facilitate the teacher's control on the student, and allowing the access to different work levels.
 This unit allows the 30 students of the classroom to visualize simultaneously all the results and the manipulation of the unit, during the process, by using a projector or an electronic whiteboard.

⑤ Cables and Accessories, for normal operation.

⑥ Manuals: This unit is supplied with 8 manuals: Required Services, Assembly and Installation, Interface and Control Software, Starting-up, Safety, Maintenance, Calibration & Practices Manuals.

Exercises and Practical Possibilities to be done with Main Items

- 1.- To use a separating calorimeter for high water contents.
- 2.- To use a throttling calorimeter for high vapour contents.
- 3.- Determination of the dryness fraction of steam.
- 4.- Global energy balance in the exchanger and study of losses.
- 5.- Determination of the exchanger efficiency. NTU method.

Additional practical possibilities:

- 6.- Sensors calibration.

Other possibilities to be done with this Unit:

- 7.- Many students view results simultaneously.
To view all results in real time in the classroom by means of a projector or an electronic whiteboard.
 - 8.- Open Control, Multicontrol and Real Time Control.
This unit allows intrinsically and/or extrinsically to change the span, gains; proportional, integral, derivate parameters; etc, in real time.
 - 9.- The Computer Control System with SCADA allows a real industrial simulation.
 - 10.- This unit is totally safe as uses mechanical, electrical and electronic, and software safety devices.
 - 11.- This unit can be used for doing applied research.
 - 12.- This unit can be used for giving training courses to Industries even to other Technical Education Institutions.
 - 13.- Control of the TCESC unit process through the control interface box without the computer.
 - 14.- Visualization of all the sensors values used in the TCESC unit process.
- By using PLC-PI additional 19 more exercises can be done.
- Several other exercises can be done and designed by the user.

a) Industrial configuration

⑦ **PLC. Industrial Control using PLC** (it includes PLC-PI Module plus PLC-SOF Control Software):

-PLC-PI. PLC Module:

Metallic box.
 Circuit diagram in the module front panel.
 Digital inputs(X) and Digital outputs (Y) block: 16 Digital inputs. 14 Digital outputs.
 Analog inputs block: 16 Analog inputs.
 Analog outputs block: 4 Analog outputs.
 Touch screen.
 Panasonic PLC:

High-speed scan of 0.32 µsec. Program capacity of 32 Ksteps. High-speed counter. Multi-point PID control.
 Digital inputs/outputs and analog inputs/outputs Panasonic modules.

-TCESC/PLC-SOF. PLC Control Software:

For this particular unit, always included with PLC supply.

Practices to be done with PLC-PI:

- 1.- Control of the TCESC unit process through the control interface box without the computer.
- 2.- Visualization of all the sensors values used in the TCESC unit process.
- 3.- Calibration of all the sensors included in the TCESC unit process.
- 4.- Hand on of all the actuators involved in the TCESC unit process.
- 5.- Realization of different experiments, in automatic way, without having in front the unit. (This experiment can be decided previously).
- 6.- Simulation of outside actions, in the cases hardware elements do not exist. (Example: test of complementary tanks, complementary industrial environment to the process to be studied, etc).
- 7.- PLC hardware general use and manipulation.
- 8.- PLC process application for TCESC unit.
- 9.- PLC structure.
- 10.- PLC inputs and outputs configuration.
- 11.- PLC configuration possibilities.
- 12.- PLC programming languages.
- 13.- PLC different programming standard languages.
- 14.- New configuration and development of new process.
- 15.- Hand on an established process.
- 16.- To visualize and see the results and to make comparisons with the TCESC unit process.
- 17.- Possibility of creating new process in relation with the TCESC unit.
- 18.- PLC Programming exercises.
- 19.- Own PLC applications in accordance with teacher and student requirements.

b) Technical and Vocational Education configuration

⑧ **TCESC/CAI. Computer Aided Instruction Software System.**

This complete package consists on an Instructor Software (INS/SOF) totally integrated with the Student Software (TCESC/SOF).

-INS/SOF. Classroom Management Software (Instructor Software):

The Instructor can:

- Organize Students by Classes and Groups.
- Create easily new entries or delete them.
- Create data bases with student information.
- Analyze results and make statistical comparisons.
- Print reports.
- Develop own examinations.
- Detect student's progress and difficulties.

-TCESC/SOF. Computer Aided Instruction Software (Student Software):

It explains how to use the unit, run the experiments and what to do at any moment.

This Software contains:

- Theory.
- Exercises.
- Guided Practices.
- Exams.

⑨ **TCESC/FSS. Faults Simulation System.**

Faults Simulation System (FSS) is a Software package that simulates several faults in any EDIBON Computer Controlled Unit.

The "FAULTS" mode consists on causing several faults in the unit normal operation. The student must find them and solve them.

There are several kinds of faults that can be grouped in the following sections:

Faults affecting the sensors measurement:

- An incorrect calibration is applied to them.
- Non-linearity.

Faults affecting the actuators:

- Actuators channels interchange at any time during the program execution.
- Response reduction of an actuator.

Faults in the controls execution:

- Inversion of the performance in ON/OFF controls.
- Reduction or increase of the calculated total response.
- The action of some controls is annulled.

On/off faults:

- Several on/off faults can be included.

c) Higher Education and/or Technical and Vocational Education configuration

⑩ **TCESC/CAL. Computer Aided Learning Software (Results Calculation and Analysis).**

This Computer Aided Learning Software (CAL) is a Windows based software, simple and very easy to use.

CAL is a class assistant that helps in making the necessary calculations to extract the right conclusions from data obtained during the experimental practices.

CAL will perform the calculations.

CAL computes the value of all the variables involved.

It allows to plot and print the results. Between the plotting options, any variable can be represented against any other.

Different plotting displays.

It has a wide range of information, such as constant values, unit conversion factors and integral and derivative tables.

d) Multipost Expansions options

⑪ **Mini ESN. EDIBON Mini Scada-Net System.**

EDIBON Mini Scada-Net System allows up to 30 students to work with a Teaching Unit in any laboratory, simultaneously.

The Mini ESN system consists on the adaptation of any EDIBON Computer Controlled Unit with SCADA integrated in a local network.

This system allows to view/control the unit remotely, from any computer integrated in the local net (in the classroom), through the main computer connected to the unit.

Main characteristics:

- It allows up to 30 students to work simultaneously with the EDIBON Computer Controlled Unit with SCADA, connected in a local net.
- Open Control + Multicontrol + Real Time Control + Multi Student Post.
- Instructor controls and explains to all students at the same time.
- Any user/student can work doing "real time" control/multicontrol and visualisation.
- Instructor can see in the computer what any user/student is doing in the unit.
- Continuous communication between the instructor and all the users/students connected.

Main advantages:

- It allows an easier and quicker understanding.
- This system allows you can save time and cost.
- Future expansions with more EDIBON Units.

The system basically will consist of:

This system is used with a Computer Controlled Unit.

- Instructor's computer.
- Students' computers.
- Local Network.
- Unit- Control Interface adaptation.
- Unit Software adaptation.
- Webcam.
- Mini ESN Software to control the whole system.
- Cables and accessories required for a normal operation.

*Specifications subject to change without previous notice, due to the convenience of improvements of the product.



C/ Del Agua, 14. Polígono Industrial San José de Valderas.
28918 LEGANÉS. (Madrid). SPAIN.

Phone: 34-91-6199363 FAX: 34-91-6198647

E-mail: edibon@edibon.com WEB site: www.edibon.com

Issue: ED01/13
Date: August/2013

REPRESENTATIVE: